

Agarose Gel DNA Loading Dye, 6x

Ultrapure Grade

Description

ROJETechnologies Agarose Gel Loading Dye is designed to optimize loading nucleic acids in agarose gels by imparting color to the sample to ensure efficient sample distribution into each well. It is provided in ready-to-use form.

specification

- Convenient ready-to-use.
- Store at 4°C.
- Dyes: Bromophenol blue (in Hashin loading dye), Orange G (in TuTik loading dye), Cresol red (in Sor loading dye).
- Due to chemical material usage that may be hazardous, always make sure to wear suitable lab coat and disposable gloves.

Protocol

Loading Dye is supplied as a 6x concentrate; thus 1 volume of Loading Dye should be added to 5 volumes of purified DNA or PCR Product.